Yan Lab by Yuntian Guan

Dobutamine Stress Echocardiography.v1

This is a protocol for assessment of cardiac function via echocardiography with dobutamine as a stress to mimic exercise. The dose of dobutamine (2.5 mg/kg) is based on many published data ranging from 1-10 mg/kg and our testing.

Procedures:

- 1. Perform baseline M-mode echo following the echocardiography protocol.
- 2. Inject the mouse with a dose of 2.5 mg/kg (10 μ l/g body weight, i.p.) dobutamine working solution (0.25 mg/ml) and start a stopwatch to track time. For example, a 30 g mouse will be injected with 300 μ l of Dobutamine working solution. Hence, 0.3 ml x 0.25 mg/ml /0.03 kg = 2.5 mg/kg.
- 3. Acquire an echo image every minute for 15 minutes (15 images post dobutamine plus one baseline image).
- 4. Follow the echocardiography protocol to export and analyze the images.

Reagent Preparation:

Dobutaimine stock solution (10 mg/ml)

- 1. Add 1 ml dH₂O to the bottle containing 10 mg dobutamine hydrochloride powder (Sigma, D0676) to final concentration of 10 mg/ml).
- 2. Mix well by pipetting.
- 3. Aliquot in 50 μ l-aliquots (500 μ g each) in light-protected Eppendorf tubes with clear labels for the date, volume, concentration and initials of the preparer.
- 4. Store the dobutamine stock solution in -70 °C freezer. Stable for at least 6 months.

Dobutamine working solution (0.25 mg/ml)

- 1. Add 1950 μ L of saline to a 50 μ L-aliquot of Dobutamine stock solution to make Dobutamine working solution (0.25 mg/ml) right before use.
- 2. Keep the Dobutamine working solution at room temperature with aluminum foil protection between the use for each mice.

References:

- Täng MS, Råmunddal T, Lindbom M, Omerovic E. Native cardiac reserve predicts survival in acute post infarction heart failure in mice. Cardiovasc Ultrasound. 2007 Dec 2;5:46. doi: 10.1186/1476-7120-5-46. PMID: 18053159; PMCID: PMC2217517.
- Heinonen SE, Merentie M, Hedman M, Mäkinen PI, Loponen E, Kholová I, Bosch F, Laakso M, Ylä-Herttuala S. Left ventricular dysfunction with reduced functional cardiac reserve in diabetic and nondiabetic LDL-receptor deficient apolipoprotein B100-only mice. Cardiovasc Diabetol. 2011 Jun 30;10:59. doi: 10.1186/1475-2840-10-59. PMID: 21718508; PMCID: PMC3141395.
- Wojewoda M, Tyrankiewicz U, Gwozdz P, Skorka T, Jablonska M, Orzylowska A, Jasinski K, Jasztal A, Przyborowski K, Kostogrys RB, Zoladz JA, Chlopicki S. Exercise capacity and cardiac hemodynamic response in female ApoE/LDLR(-/-) mice: a paradox of preserved V'O2max and exercise capacity despite coronary atherosclerosis. Sci Rep. 2016 Apr 25;6:24714. doi: 10.1038/srep24714. PMID: 27108697; PMCID: PMC4842974.
- Tong CW, Stelzer JE, Greaser ML, Powers PA, Moss RL. Acceleration of crossbridge kinetics by protein kinase A phosphorylation of cardiac myosin binding protein C modulates cardiac function. Circ Res. 2008 Oct 24;103(9):974-82. doi: 10.1161/CIRCRESAHA.108.177683. Epub 2008 Sep 18. PMID: 18802026; PMCID: PMC2867079.